

Black Sea Medium (Andersen et al. 1997)

This medium is used at the NCMA for brackish water organisms. It was designed very simply by diluting f/2 or L1 medium with distilled or de-ionized water.

To prepare, add 500 mL of f/2 or L1 medium to 500 mL of de-ionized water. Autoclave.

Andersen, R.A., Morton, S.L. and Sexton, J.P. 1997. Provasoli-Guillard National Center for Culture of Marine Phytoplankton Catalog of Strains. J. Phycol. 33 (suppl.): 1-75.